

PureCol® Collagen

Catalog Number: 5005-B

OVERVIEW

PureCol®, is approximately 3 mg/mL, pH 2, available in 100 mL volume, and produced by aseptic processing. PureCol is about 97% Type I collagen with the remainder being comprised of Type III collagen. PureCol contains a high monomer content as judged by gel permeation chromatography.

STORAGE

PureCol: The product ships on frozen gel packs with storage at 2–10° C recommended. Do not freeze. The expiration date is printed on the product label and certificate of analysis for each specific lot. The expiration date is applicable when product is handled and stored as directed.

CHARACTERIZATION

Purity: PureCol is ultrapure collagen ($\geq 99.9\%$ SDS-PAGE, ~97% Type I with remainder Type III collagen). SDS-PAGE shows the typical α , β and γ banding pattern. Gradual breakdown may occur over long periods of time thus creating atypical banding patterns.

Concentration: The concentration of PureCol collagen is approximately 3.0 mg/mL. The actual concentration is printed on the product label and certificate of analysis for each specific lot.

pH: Supplied in 0.01M HCl (pH ~2.0).

Endotoxin: <0.5 EU/mL (LAL assay)

INSTRUCTIONS FOR USE

1. Slowly add 1 part of chilled 10X PBS or 10X culture media to 8 parts of chilled collagen solution with gentle swirling.
2. Adjust pH of mixture to 7.2–7.6 using sterile 0.1 M NaOH. Monitor pH adjustment carefully (pH meter, phenol red, or pH paper).
3. Adjust final volume to a total of 10 parts with sterile water.
4. To prevent gelation, maintain temperature of mixture at 2–10° C.
5. To form gel, warm to 37° C. Allow approximately 90 to 120 minutes for gel formation.

PRECAUTIONS

This product is for R&D use only and is not intended for human or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

MANUFACTURER INFORMATION

Advanced BioMatrix, Inc.
P.O. Box 502403
San Diego, CA 92150
www.advancedbiomatrix.com
U.S. Phone: (800) 883-8220
Intl. Phone: (858) 451-2421
Fax: (510) 217-3452
support@advancedbiomatrix.com

PRODUCT WARRANTY

BioTime, Inc. and/or its subsidiaries warrants its products as set forth in the General Terms and Conditions of Sale found on ESI BIO's website at www.esibio.com/termsandconditions.

©2014 BioTime, Inc. All rights reserved. ESI BIO, ESI BIO logo and all other trademarks mentioned herein are the property of BioTime, Inc. and/or its subsidiaries or their respective owners.

For support visit www.esibio.com/support or email techsupport@esibio.com