

HyStem[®]-HP Hydrogel Kit with PEGSSDA, 7.5 mL

Catalog Number: GS315P

OVERVIEW

The HyStem[®]-HP Hydrogel Kit with PEGSSDA is composed of Heprasil[®] (thiol-modified sodium hyaluronate with thiol-modified heparin), Gelin-S[®] (thiol-modified gelatin), PEGSSDA[™] (disulfide containing polyethylene glycol diacrylate), and degassed deionized water (DG Water). Solutions of Heprasil and Gelin-S form a transparent hydrogel when mixed with PEGSSDA. All lyophilized solids are blanketed by nitrogen and under a slight vacuum.

STORAGE

Heprasil and Gelin-S: Store Heprasil and Gelin-S in original vials at -20°C for up to one year. Do not uncap the Heprasil and Gelin-S vials since both materials will crosslink in the presence of oxygen. Use a syringe to add DG Water and remove product from the vials.

PEGSSDA: Store PEGSSDA in the original vial at -20°C for up to one year. Reconstituted solutions can be stored at -20°C for ~ one month.

Note: It is recommended to reconstitute each vial in its entirety.

INSTRUCTIONS FOR USE

Heprasil, Gelin-S and PEGSSDA solutions are prepared by dissolving the lyophilized solids in the DG Water. When reconstituted, they will be in 1X phosphate buffered saline (PBS), pH ~7.4. Heprasil and Gelin-S vials contain 10 mg of material each and when reconstituted according to instruction will produce a 1% (w/v) solution. PEGSSDA vials contain 10 mg of material and when reconstituted according to instructions will produce a 2% (w/v) solution.

HyStem-HP hydrogels (3 X 2.5 mL = 7.5mL) should be prepared in the following manner:

1. Allow the Heprasil, Gelin-S, PEGSSDA, and DG Water vials to come to room temperature.
2. Under aseptic conditions and using a syringe add 1.0 mL of DG Water to the Heprasil vial. Repeat for the Gelin-S vial.
3. Place both vials horizontally on a rocker or shaker. It will take ~40 minutes for the solids to fully dissolve. Warming to not more than 37°C and/or gentle vortexing will speed dissolution. Solutions will be clear and slightly viscous.
4. Under aseptic conditions and using a syringe add 0.5 mL of DG Water to the PEGSSDA vial. Invert several times to dissolve.
5. As soon as possible, but within 2 hours of making the solutions, mix equal volumes of Heprasil and Gelin-S. To mix, pipette back and forth to mix.
6. If encapsulating cells, resuspend cell pellet in 2.0 mL of Heprasil + Gelin-S. Pipette back and forth to mix.
7. To form the hydrogel, add PEGSSDA to the Heprasil + Gelin-S mix in a 1:4 volume ratio (0.5 mL PEGSSDA to 2.0 mL (Heprasil + Gelin-S) mix by pipette).
8. If encapsulating cells, allow solution to react for 10 minutes then mix again by pipette to ensure even distribution of cells.
9. Gelation will occur within ~30 minutes.

DISSOLUTION

Dissolution of gels with cells on top and encapsulated (gel volume of 0.6 mL) in a 24 well plate. The following procedure was optimized particularly for the aforementioned gel geometry. Dissolution of gels with alternate geometry and/or volumes may require adjustments to the protocol. In general, for dissolution add at least twice the gel volume of N-Acetyl-L-Cysteine.

1. Make up the appropriate amount of 40mM N-Acetyl-L-Cysteine in 1X PBS or media and pH to 7.4.
2. Add 2 mL of 40mM N-Acetyl-L-Cysteine to the top of each gel and let sit at 37°C for 1 hour. Agitation by orbital shaking will help decrease dissolution time.

3. Confirm dissolution by pipetting solution in well up and down, observing any remaining gel. If needed, allow another 30 minutes for complete gel dissolution.
4. Remove liquid from well and place in conical centrifuge tube. If necessary, add PBS to a total of 5mL of liquid.
5. Centrifuge at 1000 RPM for 5 minutes.
6. Aspirate off liquid and process cells as desired.

Note: Each kit component has been manufactured under aseptic conditions and tested for bacteria and fungus.

PRODUCT WARRANTY

BioTime, Inc. and/or its subsidiaries warrants its products as set forth in the General Terms and Conditions of Sale found on ESI BIO's website at www.esibio.com/termsandconditions.

©2014 BioTime, Inc. All rights reserved. ESI BIO, ESI BIO logo and all other trademarks mentioned herein are the property of BioTime, Inc. and/or its subsidiaries or their respective owners.

For support visit www.esibio.com/support or email techsupport@esibio.com

